

Apo-Ident settings

When the program is started for the first time, the settings open automatically. By default, a demo profile is saved, which is used for presentations. **However, you cannot create valid test reports with the demo profile!**

Report settings

Settings > Report Settings > To create your own profile, click on the Configuration profile button on the right side of the „+“ sign

Enter the name of your pharmacy as the profile name and confirm with **<OK>**.

Another window will open asking you to enter your licence key

Note: *If you use Apo-Ident in more than one pharmacy, you need a separate licence key for each pharmacy and you have to create a separate configuration profile for each pharmacy.*

For new customers, the licence key is inserted by our sales staff at the time of delivery.

Thereafter, you will find it on the desktop as a PDF under 'Licence documents' in the 'Apo-Ident' folder or on the USB flash drive supplied.

You will need your licence key again in the following cases:

- Re-installation
- Change of computer






If you have misplaced your licence key or need support, please contact our customer service on telephone +49 351 212 496 33 or via e-mail to kundenservice@apo-ident.de.

Filename starts with > Here you can select whether the „Primary substance name“ (English) or, if available, the „Latin substance name“ should be used in the file name of the test report.

Archive location > If a profile is created, the software automatically saves the archive (test reports) on the desktop under Desktop/ApoIdent/Archiv/Profile_Name1

If a second profile is created, the software also saves the second archive under Desktop/Apo-Ident/Archiv/Profil_Name2

This ensures that several profiles are not saved in one and the same archive and that no errors occur while retrieving the archive.

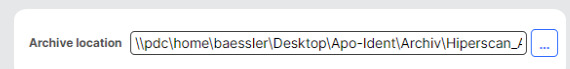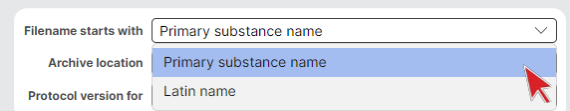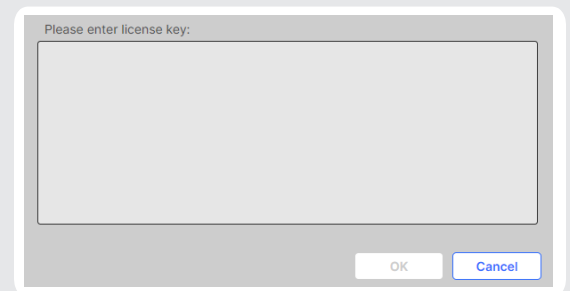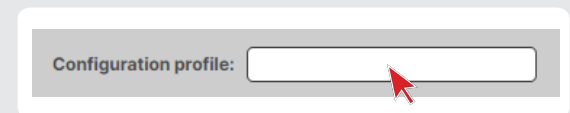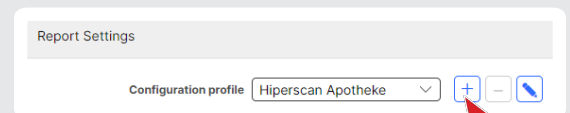

Create a configuration profile



Note: During the initial installation by our sales staff, the folder structure „Apo-Ident“ is created for you, which integrates the archive. If you would like to change the destination for saving files, first move the entire „Apo-Ident“ folder from your desktop to the new storage location. This may be a local drive or a network drive on your PC. You can change the archive directory by clicking on the folder symbol under „Profile storage location“ in Settings, Report settings. In the „Select archive directory“ window that opens, select the appropriate drive on the left and the desired folder on the right where you want to move the „Apo-Ident“ folder. Closing the settings window will transfer your changes. In the menu bar, you can use the „Archive“ button to check whether the new path has been accepted.

Report version for > The language or form of the test report for the selected profile is preset. You cannot change this.

Show difference of back projection > Section 4.3

Note: If after measuring and saving the report, you notice that the report version needs to be changed, the measuring has to be repeated after changing the necessary settings.

